# LightCycler<sup>®</sup> 480 System II

*Rapid by Nature – Accurate by Design*

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\* This product is covered in-part by US 5,871,908 or any foreign equivalents, co-exclusively licensed from Evotec OAI AG. The purchase price includes a license to practice the methods covered by US 5,871, 908 by using the product. Purchase of this product, however, does not convey to the purchaser a license or right to (i) commercially make, have made or sell reagents and/or kits, or (ii) buy or use reagents and/or kits provided by a third party used in conjunction with the product or any other thermocycler to practice the methods covered by US 5,871,908 or any foreign equivalents.

● Parts of the Software used for the LightCycler 480 System are licensed from Idaho Technology Inc., Salt Lake City, UT, USA. This product is covered by one or more of U.S. 6,197,520, 6,303,305, 6,387,621, 6,503,720, 6,730,501 and corresponding claims in their non-U.S. counterparts, owned by Roche Diagnostics GmbH and/or licensed from Idaho Technology, Inc.

▼ This LightCycler® 480 Real-Time PCR System is a real-time thermal cycler licensed for use in research under U.S. Patent No. 6,814,934 and corresponding claims in its non-U.S. counterparts, and under one or more of U.S. Patents Nos. 5,038,852, 5,656,493, 5,333,675, 5,475,610, 5,602,756, 6,703,236, 7,238,517, or corresponding claims in their non-U.S. counterparts, owned by Applera Corporation. No right is conveyed expressly, by implication, or by estoppel under any other patent claim, such as claims to apparatus, reagents, kits, or methods such as 5' nuclease methods. This instrument is for research use only. For further information on purchasing licenses for research and other non-in vitro diagnostic applications, contact the Director of Licensing at Applied Bio systems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

✦ The purchase price of this product includes a limited, nontransferable license under U.S. Patent Nos. 5,994,056 and 6,171,785 and corresponding patent claims outside the United States, owned by Roche Molecular Systems, Inc. and F. Hoffmann-La Roche Ltd ( "Roche" ), and under U.S. Patent No. 6,569,627 and corresponding patent claims outside the United States, licensed from Idaho Technology Inc., for using only this amount of the product for dsDNA-binding dye processes covered by said patent solely for the purchaser's own internal research and development activities. This product is also a Licensed Dye Binding Kit for use with service sublicenses available from Applied Biosystems. No right under any other patent claims (such as apparatus or system claims in U.S. Patent No. 6,814,934) and no right to use this product for any other purpose or for commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is hereby granted expressly, by implication, or by estoppel. This

product is for research use only. Diagnostic uses require a separate license from Roche.

Further information on purchasing licenses to practice real-time PCR processes may be obtained by contacting the Director of Licensing at Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

◆ A license to perform the patented 5' Nuclease Process for research is obtained by the purchase of (i) both Authorized 5' Nuclease Core Kit and Licensed Probe, (ii) a Licensed 5' Nuclease Kit, or (iii) license rights from Applied Biosystems. This product is an Authorized 5' Nuclease Core Kit. Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 5,079,352, 5,789,224, 5,618,711, 6,127,155, 5,677,152, 5,773,258, 5,407,800, 5,322,770, 5,310,652, 5,210,015, 5,487,972, and claims outside the US corresponding to US Patent No. 4,889,818. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. Separate purchase of a Licensed Probe would convey rights under the applicable claims of US Patents Nos. 5,538,848, 5,723,591, 5,876,930, 6,030,787, 6,258,569, 5,804,375 (claims 1-12 only), and 6,214,979, and corresponding claims outside the United States. No right under any other patent claim and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained from the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

✦ The purchase price of this product includes a limited, nontransferable license under U.S. Patent Nos. 6,174,670, 6,245,514 and corresponding patent claims outside the United States, licensed from Idaho Technology Inc., to use only this amount of the product for HybProbe assays and related processes described in said patents solely for the research and development activities of the purchaser.

No right under any other patent claims (such as apparatus or system claims in U.S. Patent No. 6,814,934) and no right to use this product for any other purpose or for commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is hereby granted expressly, by implication or estoppel.

This product is for research use only. Diagnostic uses require a separate license from Roche.

Further information on purchasing licenses to practice real-time PCR processes may be obtained by contacting the Director of

Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

❖ A license to perform the patented 5' Nuclease Process for research is obtained by the purchase of (i) both Licensed Probe and Authorized 5' Nuclease Core Kit, (ii) a Licensed 5' Nuclease Kit, or (iii) license rights from Applied Biosystems. This product contains Licensed Probe. Its purchase price includes a limited, non-transferable immunity from suit under U.S. Patents Nos. 6,214,979 and 5,804,375 (claims 1-12 only) and corresponding patent claims outside the United States. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. The right to use this product in the 5' Nuclease Process under the applicable claims of US Patents Nos. 5,210,015 and 5,487,972, and corresponding patent claims outside the United States, can be obtained through purchase of an Authorized 5' Nuclease Core Kit. Except under separate license rights available from Applied Biosystems, no right under any other patent claim, or to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, or to sublicense, repackage with other products, or resell in any form, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained from the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

○ The technology used for the LightCycler<sup>®</sup> System is licensed from Idaho Technology Inc., Salt Lake City, UT, USA.

● ProbeLibrary is covered by US and other patent applications owned by Exiqon A/S. Locked Nucleic Acids (LNA) are covered by U.S. Patents No US 6,794,499, US 6,670,461, US 6,268,490 & US 6,770,748 and other patents and patent applications owned by Exiqon A/S and Prof. Takeshi Imanishi. The quencher used in the probes is covered by patent applications owned by Exiqon A/S.

★ This product is sold for use by purchaser only. The sale or distribution of this product or any product derived from the use of this product to any third party is expressly excluded. Especially excluded is the commercial use or sale of oligonucleotides or other products labeled with this product.

+ Under license from Gilead Sciences Inc.

◀ No license is conveyed with the purchase of this product under any of US Patents Nos. 5,210,015, 5,487,972, 5,804,375, 5,994,056, 6,171,785, 6,214,979, 5,538,848, 5,723,591, 5,876,930, 6,030,787, and 6,258,569, and corresponding patents outside the United States, or any other patents or patent applications, relating to the 5' Nuclease and dsDNA-Binding Dye Processes. For further information contact the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

◇ Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 5,079,352, 5,789,224, 5,618,711, 6,127,155, 5,677,152, 5,773,258, 5,407,800, 5,322,770, 5,310,652, 5,994,056, 6,171,785, and claims outside the US corresponding to US Patent No. 4,889,818. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim (such as apparatus or system claims in US Patent No. 6,814,934) and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

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LIGHTCYCLER, HYBPROBE, and SIMPLEPROBE are trademarks of Roche. SYBR is a registered trademark of Molecular Probes, Inc. PROBELIBRARY and LNA are registered trademarks of Exiqon A/S, Vedbaek, Denmark. Other brands and product names are trademarks of their respective holders.

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